


How To Peak evaluation with Chip DSC

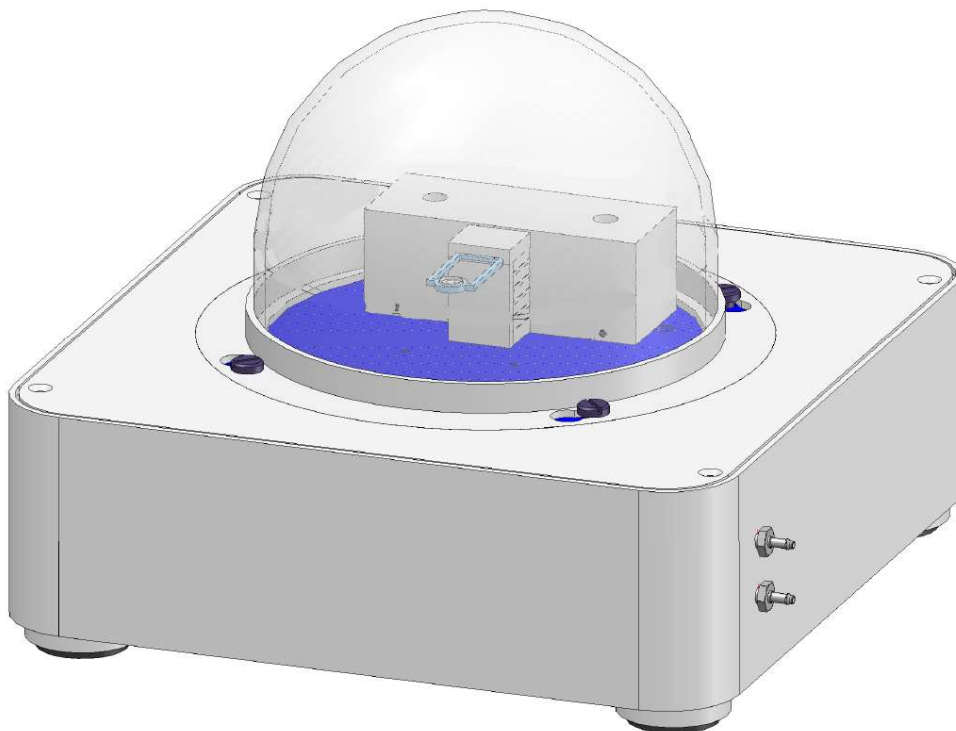

Linseis Messgeräte GmbH

Gerlach

Date: 13.10.2020

Index

1. General information.....	3
2. Manual peak evaluation.....	3
3. Double peak evaluation	4
4. Automatic peak evaluation	5

1. General information


This manual is a short description of how to evaluate peaks of Chip-DSC measurements. The evaluation of peaks, beginning with the type of used baseline up to setting initial and endpoint is important to obtain valid results. To receive reproducible values clear peaks and measurements without distortion are required. For more Information, read the other available instructions about software or specific manuals for the Chip DSC.

2. Manual peak evaluation

- Load and select curve
- Choose peak evaluation tool




- Choose baseline character

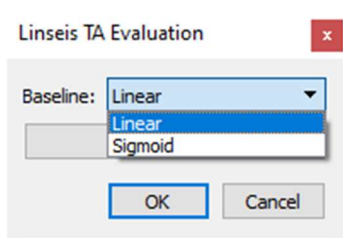


Linear

linear baseline between two points

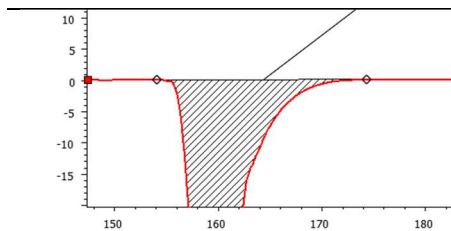


Sigmoid

baseline by two baselines before and after transition

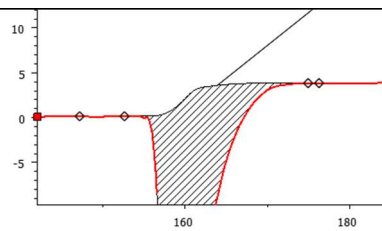

- Select information to your needs

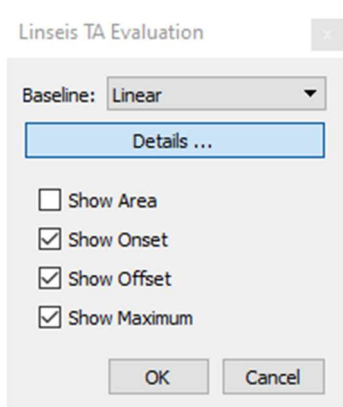


3. Double peak evaluation


- Load and select curve
- Choose peak evaluation tool



- Evaluate first peak

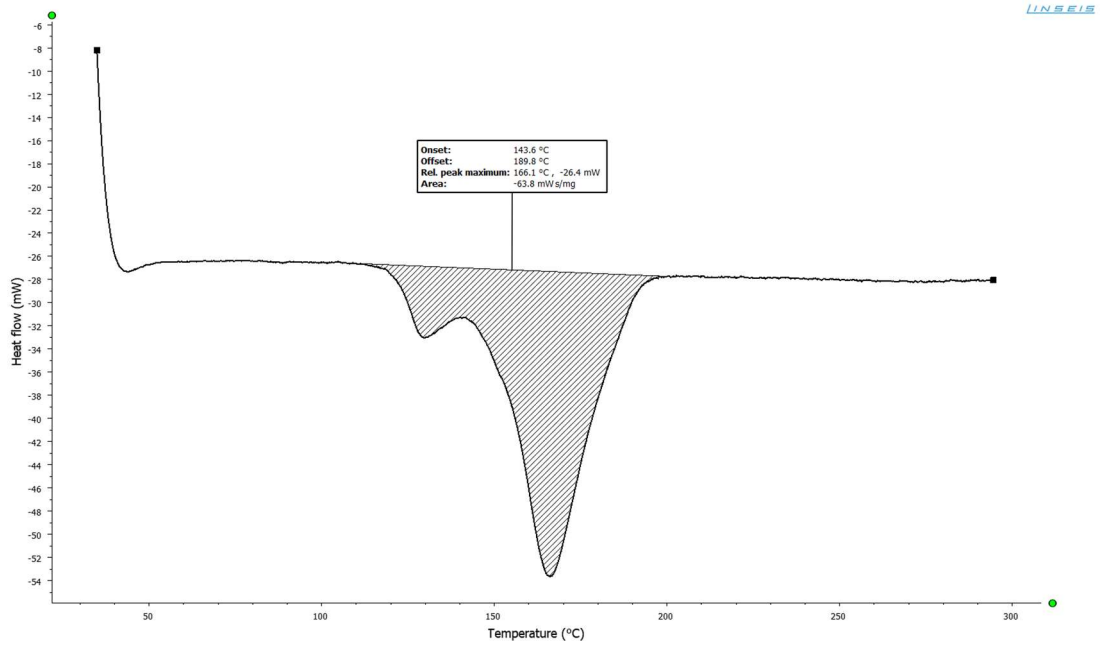

- Evaluate second peak

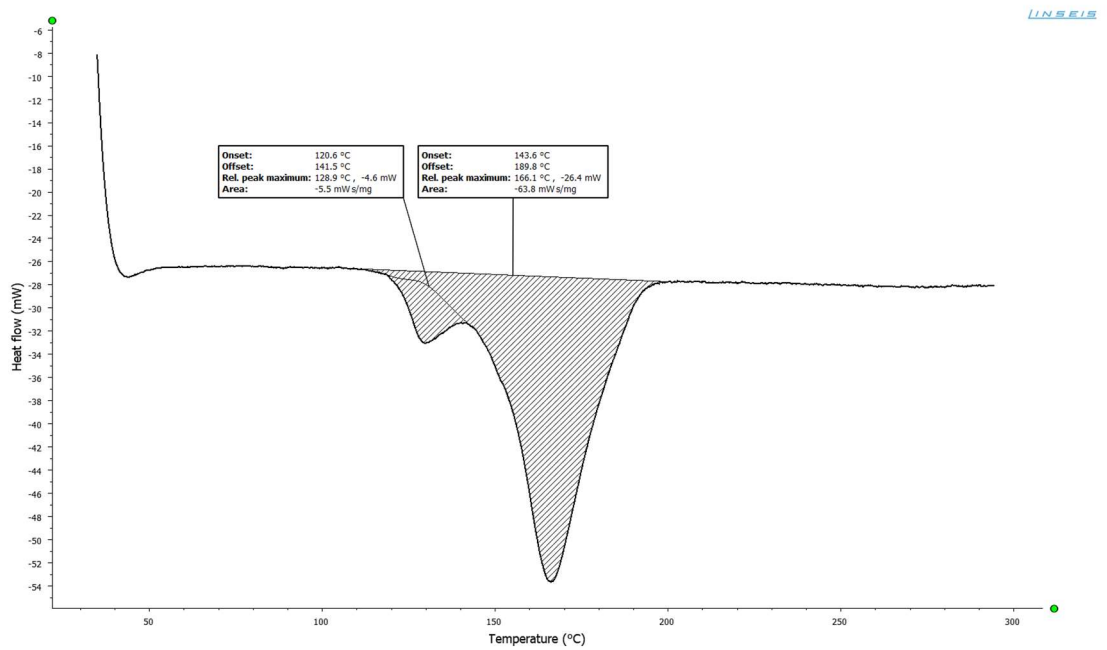


- Subtract peaks

4. Automatic peak evaluation

- Load and select curve
- Choose automatic peak evaluation tool



- Check received information

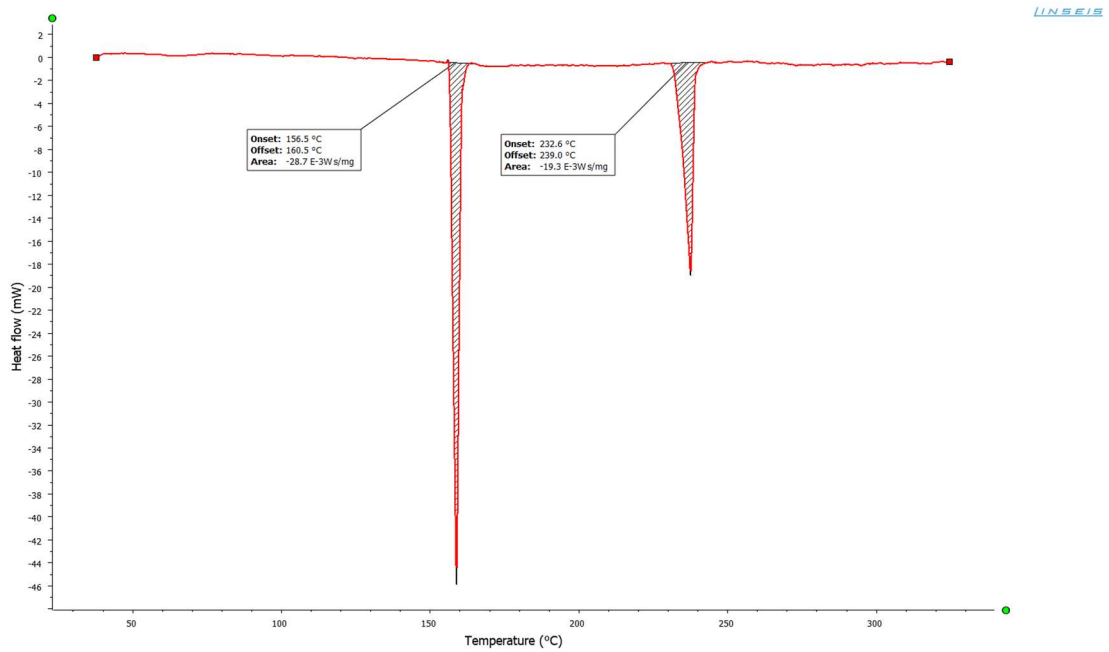

- Adjust peak settings (double click on label)

