

# How To Setup your Linseis PC



Linseis Messgeräte GmbH  
Gerlach  
Stand: 10.07.2019

## Index

1. Setup the hardware.....	3
2. Setup Windows OS .....	3
3. Setup the Linseis Software .....	5

## 1. Setup the hardware

- At first make sure everything is delivered you need to setup your PC correctly



<input type="checkbox"/>	Tower
<input type="checkbox"/>	Monitor
<input type="checkbox"/>	Keyboard
<input type="checkbox"/>	Mouse
<input type="checkbox"/>	Power cables
<input type="checkbox"/>	Video cable
<input type="checkbox"/>	USB-Stick (you can find this with your measuring device)

- Connect your monitor with the tower
- Connect mouse and keyboard with the tower
- Connect monitor and tower to power supply and turn everything on

## 2. Setup Windows OS

- Turn the PC on



- For the first time Windows will setup itself, you just need to follow the instructions and insert the information to your needs


- You need to connect the PC with the internet while setting up the device (Ethernet or WIFI)



- When the WINDOWS installation is finished, connect the provided USB-Stick with the tower




- Start the "Install"-Batch file and follow the instructions in the CMD-Window

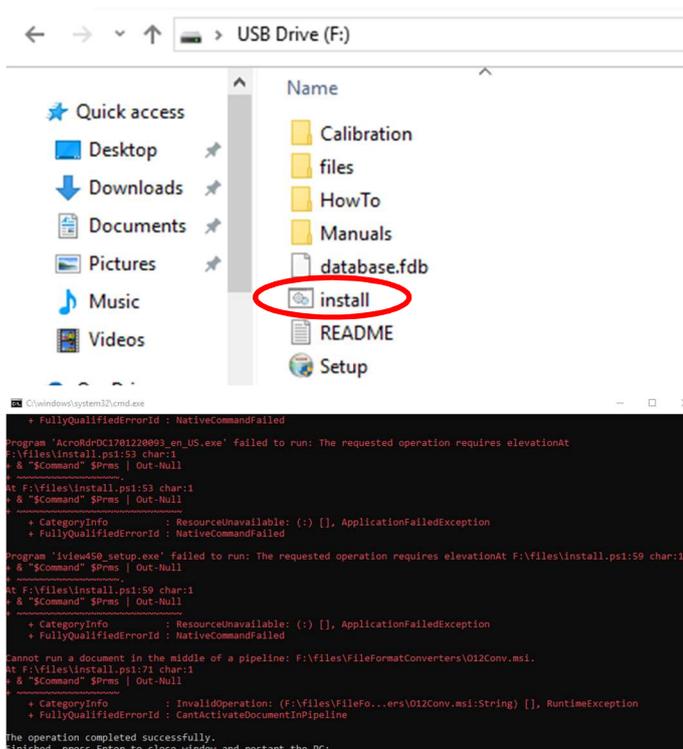


- After you restart the PC, everything is setted up on your brand new Linseis PC

## 3. Setup the Linseis Software

- If the Windows-Setup finished you can start install the Linseis Software on your PC
- Start the Setup.exe from your USB-Stick

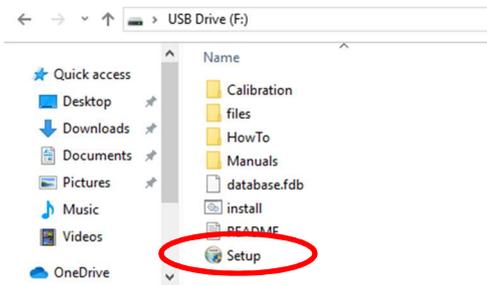


- Choose the device you want to install on your PC in the dropdown menu

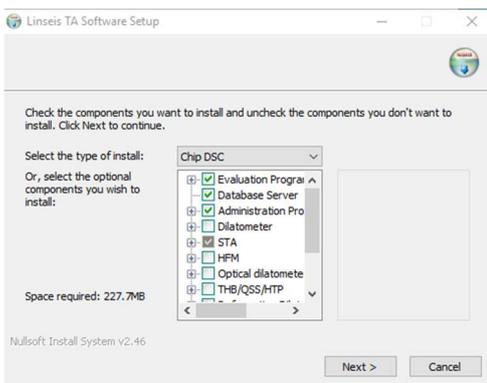

- Click Ok and confirm the upcoming questions

The database for your device will be copied automatically from the USB-Stick

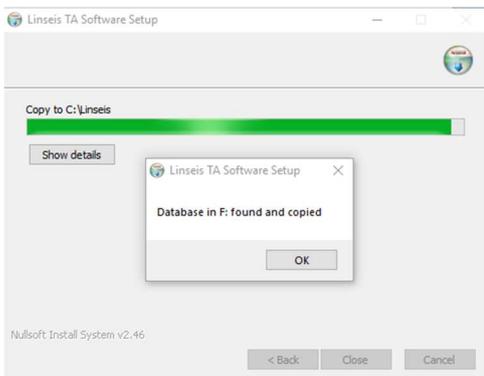

- Connect your Linseis device with the PC and follow the instructions in the manual
- Now you can Use your Linseis Device and PC to your needs

